

# specbos 1211 Firmware Changelog

## V3.4.12

- new parameter to signalize, that straylight correction matrix is to be used

## V3.4.11

- bug in new fault pixel function -> fixed

## V3.4.10

- neighboring fault pixel were not possible -> fixed

## V3.4.9

- reference measurements with integration time at split time were not correct -> fixed

## V3.4.8

- \*meas:compdark now uses the average setting instead of fixed average=1

## V3.4.7

- unexpected behavior, when pulselength of a flash light was set to more than 2200us (\*para:flash) -> fixed

## V3.4.6

- Sync measurement was not using integration times down to min\_tint -> fixed

## V3.4.4

- Sync measurement was not working (since V3.4.0) -> fixed

## V3.4.0

- Support for new detector S11155