

Updating the Firmware of JETI SPECFIRM-based devices

This documentation is valid for the following devices and electronics:


- spectraval 1501/1511
- specbos 2501
- SDCM3/4
- PE60_2
- PE65

Please note that in rare cases a firmware update may fail and the device or the electronics may become unusable.


The update should only be carried out in justified cases and is at your own risk.

Instructions for the update



1. Download the corresponding update package and unpack it.
2. Start the programming tool SFProg.exe

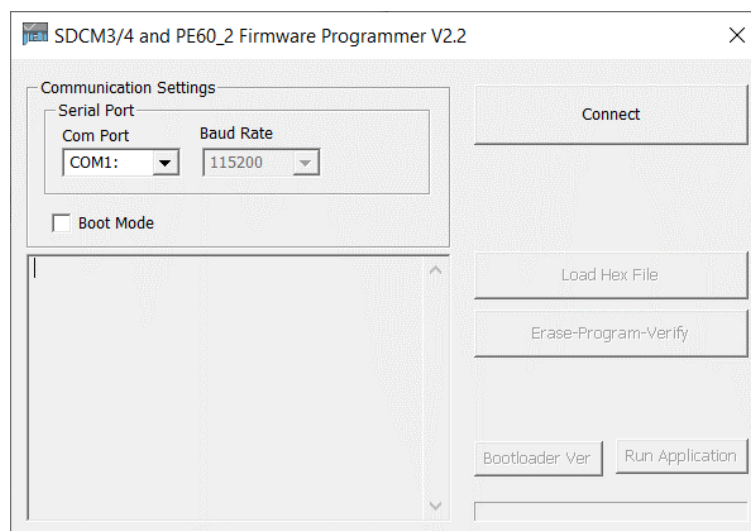


3. Connect the JETI unit via USB port to the PC
4. Select the correct COM port (in case of problems: check the COM port number in the device manager)


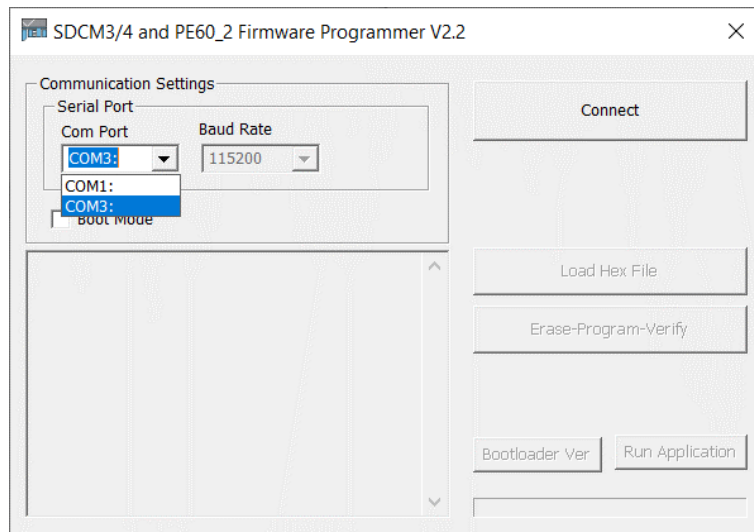


5. Load the firmware file by pressing the **Load Hex File** button and selecting it in the appropriate folder

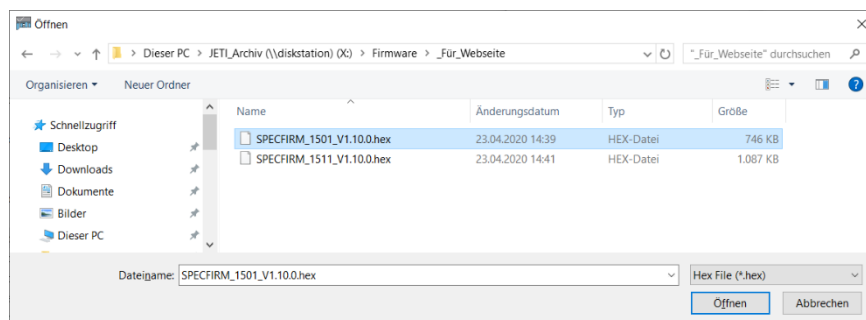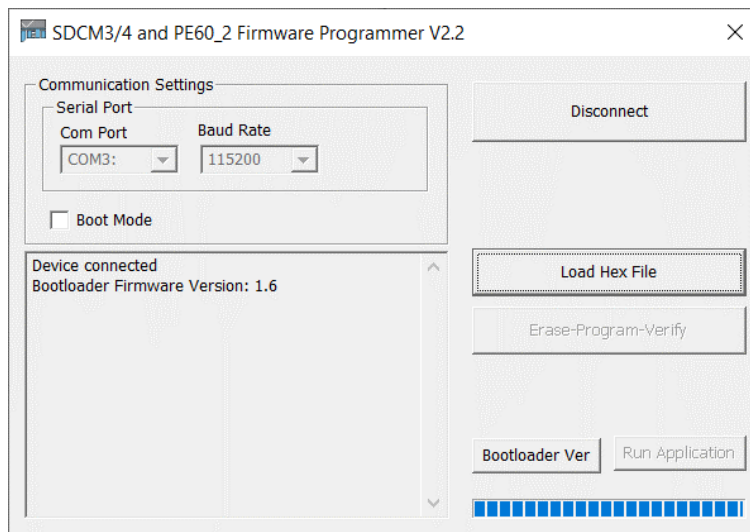

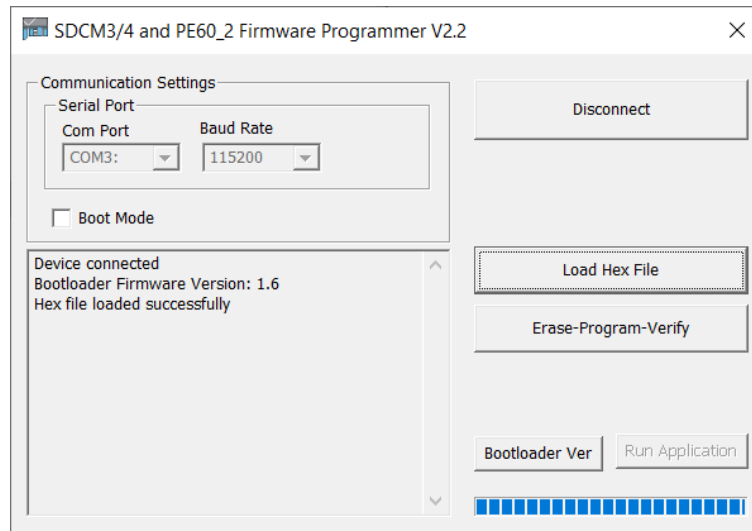

6. Press the Erase-Program-Verify button and wait for the finalization of the programming (Verification successful)

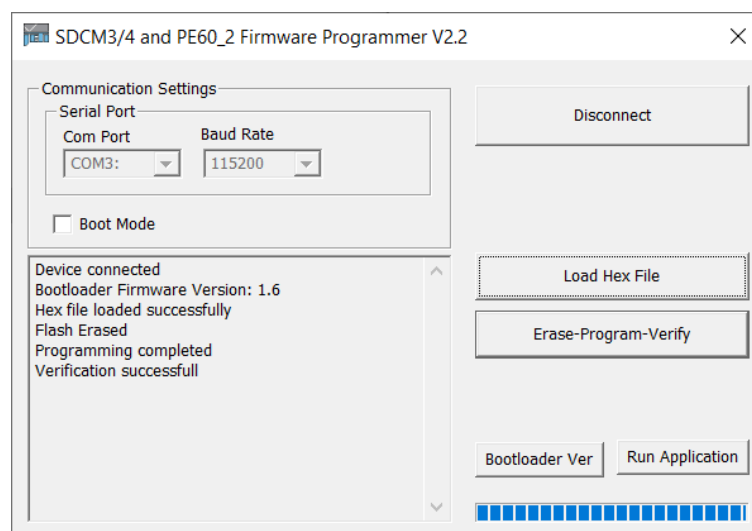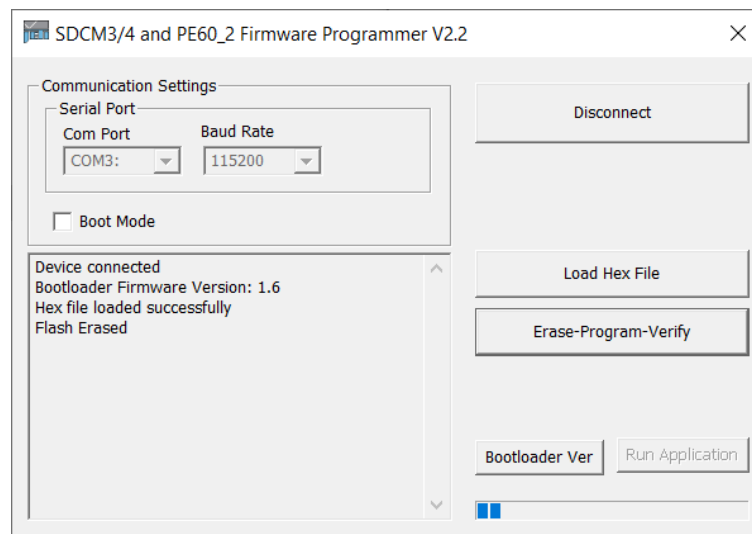

7. Press the Disconnect button, close the software and confirm the exit

